



Ingredients Notes on safety and use

1. General usage information:

Use the device only in the manner described in the manual. Do not modify the components, it results in the termination of the warranty. This could lead to a malfunction. In turn they can lead to e.g. electrocution electric current. The manufacturer is not responsible for any damage to health resulting from improper use of the equipment, reagents or any other components.

The device is not suitable for use by children and adolescents.

The durability of the reagents is confirmed by the manufacturer for one year. Subsequent use causes poor test results.

2. Reagents and other chemical equipment:

Under no circumstances should you leave open vials, bottles with reagents or powders. Pictograms for the use of chemicals:



3. Detailed information on the use of chemical reagents, methods of storage, neutralization:

a) Store reagents in:

- **Closed packages**
- In the original collective packaging



- Out of the reach of children, light, water, high temperature

b) When using the reagents:

- Keep a distance of at least several centimeters from the nose, mouth, eyes - Do not inhale the vapors

- Be careful with the skin, some substances can cause an allergic reaction

- **Do not eat !!! - if swallowed, immediately contact a doctor or the nearest poison department. Contact to the manufacturer: 699713471, orders@diodeprobe.com**

- Watch out for eyes, open wounds - in the event of the substance getting into eyes, rinse immediately and contact a doctor or the nearest poison department

- In the event of the product getting on the skin, watch for an allergic reaction, erythema, redness. In this case, contact your doctor.

- Clothing - pay special attention to care for it. Substances in solutions and powders do not damage clothing, but over time may dry out and completely destroy the material. You should rinse the clothes at least several times with water. In the case of colored compounds, after washing with water, it is recommended to wash with water with washing powder / soap.

- Neutralization of spilled preparation or scattered powder: remove substances by first dry wiping and then wet wiping several times. Rinse the surface with water.

- Disposal of old reagents, reagents expire after 1 year: remove the compound to the sewage system and rinse with a large volume of water or take it to a special collection point for used chemicals and medicines.



4. Responsibility of the manufacturer:

The manufacturer is not responsible: for any health detriment resulting from improper use of reagents, for the above-mentioned consequences of accidents. **Please store and use the reagents carefully !!!**

